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www.steinbrenner-laborsysteme.de -

Operating Manual Liquidator⁹⁶ RT

Version Dec, 2008, subject to alterations

Dear Customer,

Liquidator⁹⁶ is an easy-to-handle, fully manually operated pipetting system. 96 samples can be handled at a time making Liquidator⁹⁶ an ideal tool for high throughput applications.

IMPORTANT *Liquidator⁹⁶ is intended for use by persons knowledgeable in safe laboratory practices. Please operate the instrument according to this manual in order to guarantee safe and undisturbed long term use and display the manual close to the device. Make sure that all general and work-specific safety instructions are being followed closely.*

I. OPERATIONAL USE

Liquidator⁹⁶ is an air replacement pipettor with 96 separate channels and the 96 tip cones arranged in an 8 by 12 pattern compatible with the international microtitre plate standard. The pipetting head can be moved along a vertical and a horizontal axis and there are two movable trays each with two working positions in SBS format. Liquidator⁹⁶ works with disposable pipette tips and covers a volume range from 1 to 200 µl using the two tip sizes available. The working position in the back of the right tray is designed to hold the tip box, but has an SBS footprint also. This means that a maximum of four different plates with SBS footprint can be held simultaneously if the tip box is replaced. The functionality of the device is very similar to that of a handheld pipettor which makes switching to Liquidator an easy step for experienced labworkers.


- **Liquidator is dedicated for research purposes only.** Steinbrenner Laborsysteme GmbH will not be liable for any harm or damage resulting from using it for other purposes without their written consent.

II. PREPARATIONS FOR USE

Liquidator⁹⁶ should be placed level in a dry and clean place without exposure to sunlight. The recommended operation temperature is 20° C - 25° C. Exceeding this range may cause malfunction of the device and will lead to a loss of precision, as well as sudden changes in temperature or humidity.

- **Before starting to operate the device, make sure that the safety lock of the pipetting head is released.** Pull the black knob of the security bolt (5) on the right side of the device and turn it about 10-15 degrees such that it locks in pulled position.
- **The pipetting head has to be secure locked again prior to any further moving of the device** to avoid uncontrolled movement of the pipetting head during transportation. Move the pipetting head to the right and downwards to the end position. Keep it in this position while pulling the black knob of the security bolt (5) on the right side of the device and turn it about 15 degrees such that it comes close to the housing of the device. You will hear the security bolt snap, indicating that the pipetting head is safely locked.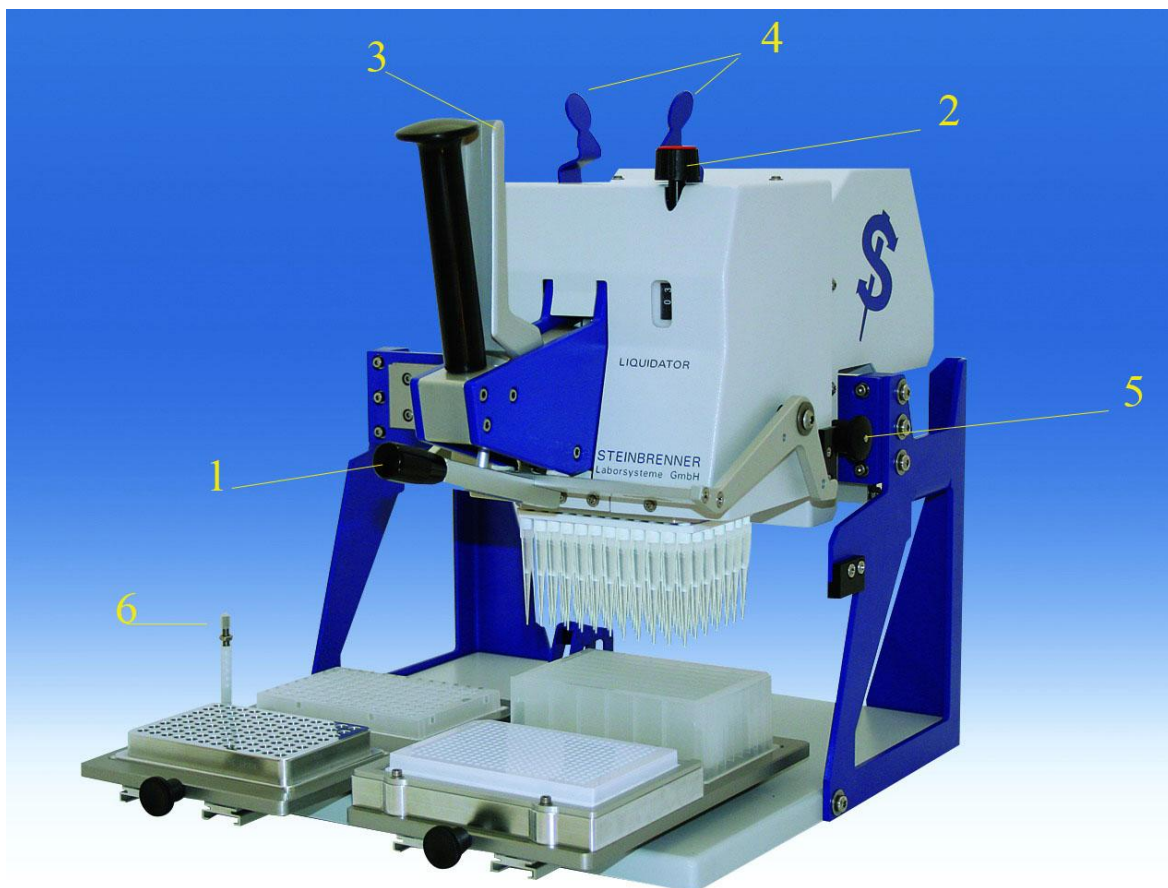

- 1 tip mounting lever
- 2 volume selector
- 3 pipetting lever
- 4 tip ejector/tip indicator
- 5 security lock
- 6 height adjuster

III. PIPETTING


Liquidator⁹⁶ works similar to conventional manual pipettors and can be operated just as easily. It requires only a few steps to efficiently work with the device. Special emphasis has been placed on an easy and comfortable handling of Liquidator⁹⁶.

Liquidator is designed for convenient and relaxed operation. Please contact us immediately if any function of your Liquidator⁹⁶ should require more force than usual.

1. Mounting of Tips

- a) Move the pipetting head smoothly upwards and to its left end position. Pull right tray forward to its end position and insert a tip box with fresh tips into the working position of the right rear base plate (left). Press down the tip box to ensure that it is correctly positioned and secured by the positioning means of the right tray.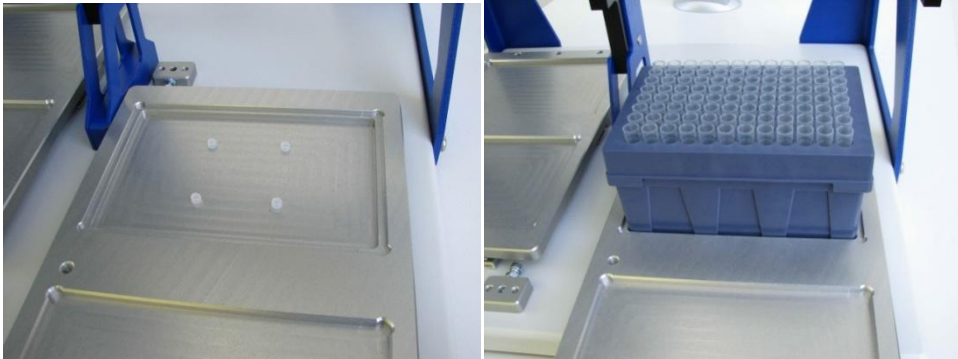

- b) Smoothly move the pipetting head to its right end position and then downward. The two grips of the tip ejection mechanism (4) move outwards, indicating that the tip cones of the pipetting head have been correctly inserted into the tips.
- c) The pipetting head can now be pulled into the tips by gently pressing downwards lever (1). *If much force is needed please check first that the tip box is correctly placed and the movable tray as well as the pipetting head are in their end positions.*

WATCH THE GRIPS (4) OF THE TIP EJECTING MECHANISM ON TOP OF THE DEVICE DURING THE PROCEDURE OF MOUNTING FRESH TIPS!

*These grips serve as a **Tip Indicator**: They move outwards when the tip cones correctly hit into the wide ends of the pipet tips of the tip box.*

- **Do not press down the tip mounting lever (1) unless the grips (4) have moved outwards a bit** when you move down the pipetting head onto the tip rack. *Sometimes it may be necessary to move the pipetting head a little bit to insert the nozzles into the tips.*

- **Stop further pressing down the tip mounting lever (1) as soon as the grips (4) have moved outwards completely.**

2. Volume setting

The standard 200 µl tips provide very good pipetting results from 3 µl to 200 µl. The volume is set and indicated by volume selector (2) while pulling pipetting lever (3) at the same time.

- **Do not screw the volume selector (2) without keeping the pipetting lever (3) pulled at the same time.**
- **Never screw volume selector (2) over 200 µl or down to less than 0 µl.**

3. Aspiration of Liquid

Gently pull the pipetting lever (3) until the first stop and keep it in this position whilst move the pipetting head down until the tips dive into the liquid. Let lever (3) slowly move back to aspirate liquid into the tips.

4. Dispensing of Liquid

- Position pipetting head over the target plate and move it down carefully. The liquid is dispensed by pulling the pipetting lever (3) to the first stop. (Multiple pulling and releasing leads to mixing of liquids and resuspension of a pellet). Liquid remaining in the pipet tips can be blown out by quickly pulling lever (3) further to its final stop.

TIPP *The vertical position of the tips can be defined by **adjustable stops** (6) individually for three of the working positions. With this feature, the distal tip end can be stopped in a vertical position without the tip ends touching the bottom of the plate. This prevents selfsealing or possible destruction of coatings or cell layers. **Please note that the adjustable means may not be used for the tip box position to avoid malfunction of the tip mounting mechanism.***

5. Tip change

- Move the pipetting head over an empty tip box and downwards until the distal ends of the tips fit into the holes of the box. The tips are ejected by pressing together the two levers (4) on top of the pipetting head.

TIPP *If you want to reuse your empty racks do not eject your used tips into a clean tip box but into another suitable container, e.g. a used deep well plate.*

IV. GENERAL PIPETTING RECOMMENDATIONS

The ideal speed of pipetting has to be found out empirically and is depending on the circumstances and the liquids to be pipetted. With the following suggestions we want to help you to facilitate this process.

PIPETTING OF SMALL AMOUNTS OF LIQUID

- In most cases the pre-rinsing of the tips - by aspirating and dispensing 2-3 times - leads to better results.
- Do not to use the blow out mode for dispensing when pipetting foaming liquids or small volumes (<10 µl)
- When pipetting small volumes it is of advantage to do this very carefully and with low speed. Touching the surface of the receiving vessel or surface of the liquid in the vessels with the ends of the tips whilst gently pipetting out may improve the result.
- Alternatively, dispense the liquid above the destination plate such that it forms a hanging drop at the distal end of the tip. By slowly approaching the bottom of a plate or the liquid surface in a plate these drops will be drawn from the tips.
- Well experienced users of Liquidator may also use the so called "tip touch"- technique: Move down the pipetting head to the height desired and *carefully dispense the liquid. Now move the tray with the destination plate until the distal ends of the tips hit the side walls of the wells.* The liquid is now drawn from the tip by capillary forces.

The „tip touch“ procedure has to be carried out very carefully to not damage the tip cones.

STERILE PIPETTING

- Liquidator⁹⁶ may also be operated in a laminar flow hood. To sterilize the surface of the device, wipe with alcohol or BARRYCIDAL, a non-toxic and non-corrosive disinfectant available from Steinbrenner Laborsysteme GmbH.

- **Expose the device to UV-radiation only as long as necessary, but never overnight!**
The radiation and the ozone generated will affect the material of the sealings and other parts.
- **Do not autoclave Liquidator⁹⁶ or parts of it and never use autoclaved tips!**
The tips as well as the boxes shrink during autoclaving, which may lead to a damage of the device! For sterile pipetting radiation-sterilized tips from Steinbrenner Laborsysteme GmbH are available.

V. MAINTENANCE AND SAFETY INSTRUCTIONS

Liquidator⁹⁶ does not require much service when used following the instructions closely. The surface can be cleaned with a soft cloth slightly moistened with water or a mild detergent. This instruction manual may not predict any situation with potential risk occurring during the use of Liquidator⁹⁶. Dangerous or harmful liquids may only be handled with suitable protective means such as safety glasses, gloves, or suits. Contaminations of the device have to be removed immediately and it fully lies in the responsibility of the user himself to follow this instruction manual and to work in accordance with the general safety guidelines set up for his institution or working area. To prevent damage and harm of personell, limits in the use of Liquidator⁹⁶ may have to be installed also.

VI. GENERAL WARNINGS

- Pipetting strong acids, bases or other aggressive liquids may damage seals, nozzles or other parts of Liquidator⁹⁶. The contact of parts of Liquidator⁹⁶ with such substances is strictly to avoid as well as the seeping of liquid into the tip cones. Accidentally applied aggressive liquids or harmful substances have to be removed immediately! Fuming or explosive liquids may not be pipetted at all.

- Liquidator⁹⁶ has to be inspected at owner's expense by Steinbrenner Laborsysteme GmbH every 2 years.
- The device is only accepted in the shipping container supplied by Steinbrenner Laborsysteme GmbH
- Liquidator⁹⁶ has to be thoroughly cleaned and disinfected/decontaminated prior to shipping with a declaration of the sender that the instrument is free of toxic or otherwise harmful substances.
- Do not work directly under the pipetting head. Move it over to have free access to the working place required.

VII. NOTICE OF DEFECTS, CLAIMS, PERIOD OF LIMITATION

Buyer undertakes to inspect the device immediately upon receipt and shall give written notice of any defects. Failure to notify Steinbrenner Laborsysteme GmbH about defects may result in Buyer's loss of right to have the defect remedied. In case of justified complaints we will remedy the defect at our discretion by repair or replacement. The period of limitation of claims arising from defects is one year from the date of delivery. Warranty claims will not be considered for defects occurring after passing of risk due to unsuitable or improper use, wrong assembly or start-up by Buyer or third parties, invalid mode of operation, normal wear and tear and improper maintenance. If Buyer or third parties rectify defects with improper remedies, Steinbrenner Laborsysteme GmbH shall not be liable for the resulting consequences. The same shall apply for any changes to the delivery item without our prior written consent.

The warranty does not extend to damages that are not product-inherent defects defined by the law or to product failure resulting from product misuse, improper surrounding conditions, or use with improper chemicals. The warranty will also be deemed as void if seals of the housing are broken or a fault is found to have been caused by e.g. maltreatment, accidental damage or use for operations outside its specified limitations as well as by using accessories such as tips, spare parts, lubricants and o-rings unauthorized by Steinbrenner Laborsysteme GmbH.

The warranty only comprises the material value as well as labour expenses to restore the function on behalf of Steinbrenner Laborsysteme GmbH. We explicitly exclude liability for financial loss, material damage and personal injuries resulting from using Liquidator⁹⁶ or resulting from a malfunction of Liquidator⁹⁶.

Liquidator⁹⁶ stays in full ownership of Steinbrenner Laborsysteme GmbH until fully paid and is sold under our Terms and Conditions of Sale which can be studied on our webpage. Alternatively, please request for a copy.

STEINBRENNER LABORSYSTEME GMBH

IN DER AU 17

D-69257 WIESENBACH

TEL. 06223-861247

FAX: 06223-861248

mail@steinbrenner-laborsysteme.de

www.steinbrenner-laborsysteme.de -