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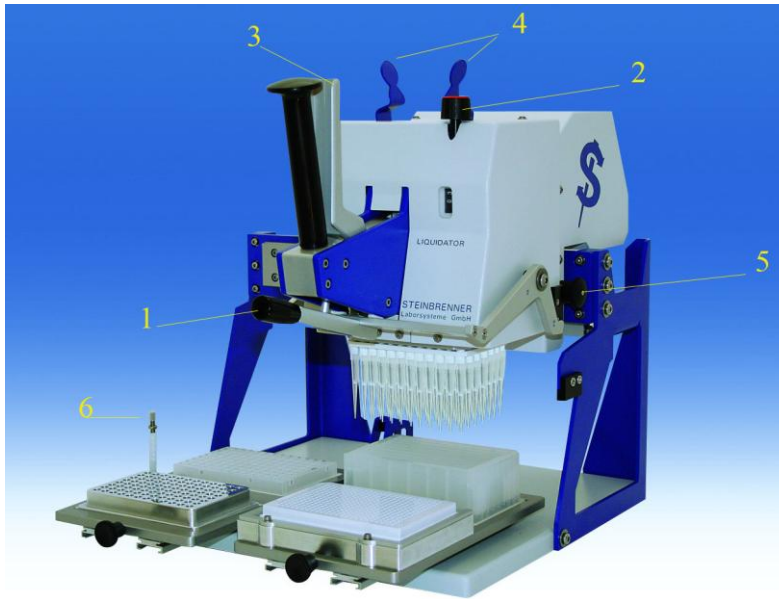
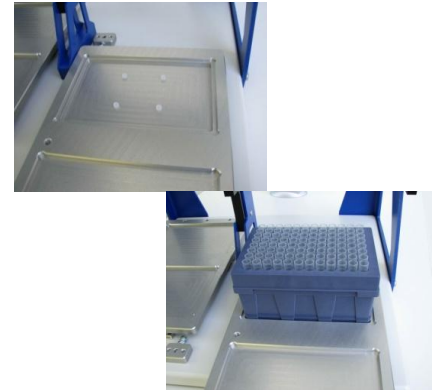
### Mounting of Tips

1. Pull right tray to the front end position. Insert tip box in the backward position and press it down until it is secured by the positioning means (below).

2. Move the pipetting head to the right end position and move it smoothly downwards over the tip rack. Watch the two tip ejector grips (4) on top of the device. By moving outwards, these grips **indicate** that the tip cones have correctly hit into the tips.

3. Press down the tip mounting lever (1)

- The tip mounting lever (1) may only be used when indicator grips (4) have moved outwards when lowering the pipetting head down onto the tip rack.
- Stop pressing downward tip mounting lever (1) as soon as you feel a click.

### Adjustment of pipetting volume

- Always keep pipetting lever (3) pulled whilst adjusting pipetting volume with screw (2).
- Never screw volume selector (2) over 200  $\mu$ l or down to less than 1  $\mu$ l.

### General Recommendations

- Liquidator is designed for convenient and relaxed operation. Please contact us immediately if any function of your Liquidator<sup>96</sup> should require more force than usual.
- Always secure the pipetting head with bolt (5) prior to moving the device to another place to prevent uncontrolled movement of the pipetting head.
- Do not autoclave Liquidator<sup>96</sup> or parts of it and never use autoclaved tips!