

Packaging Instruction


Liquidator96




Dear Customer,

Liquidator may only be shipped in the shipping box provided by Steinbrenner Laborsysteme GmbH. Please follow these instructions closely to avoid damage during transportation. Contact us prior to any shipment if you have no such shipping box or not all components required.

- 1 black foam piece (1)
- 1 cover (blue or clear) (2)
- 2 protector small (3)
- 1 protector long (4)
- 2 cable binder connected (5)
- 1 shipping box (6)
- 1 lower foam piece (7)
- 1 upper foam piece (8)
- 1 plastic bag (9)
- Safety Declaration

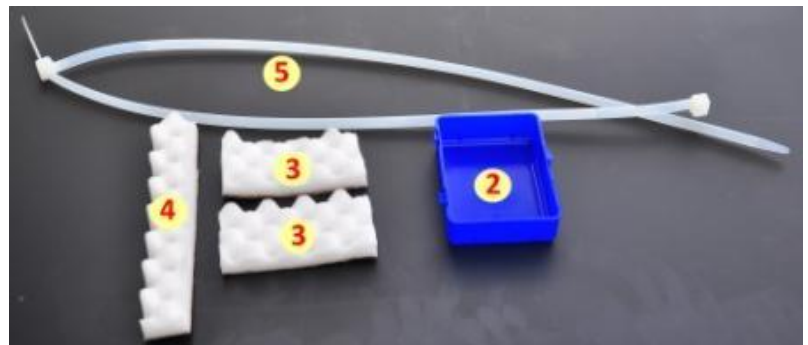


1. Move pipetting head to the centre and push cover (2) over the pin plate and hold it in this position.

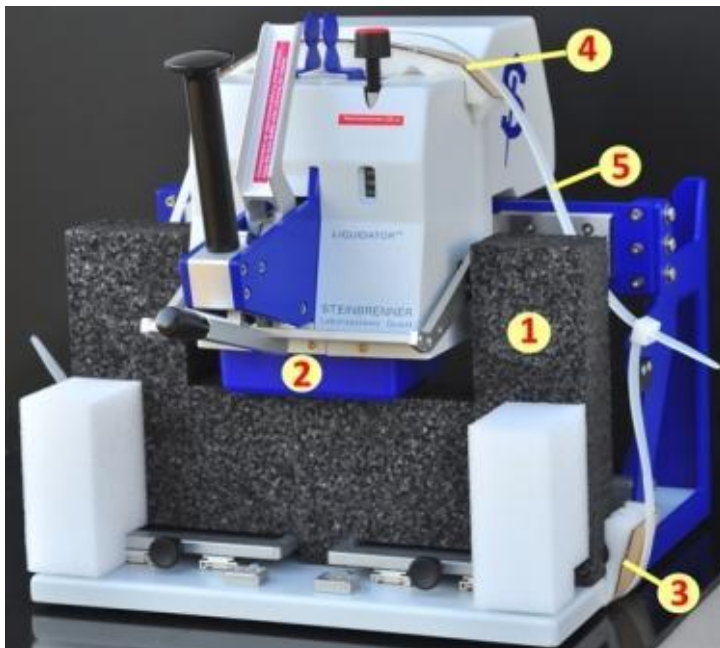

2. Insert black foam piece (1) from the front and push it between the two drawers until it rests under the pipetting head as shown, now supporting cover (2).

3. Pull cable binder around Liquidator as shown. Place protectors (3) and (4) to prevent damages of the surface of base plate and pipetting head, respectively. Make sure that the protectors do not move while pulling cable binder tight.

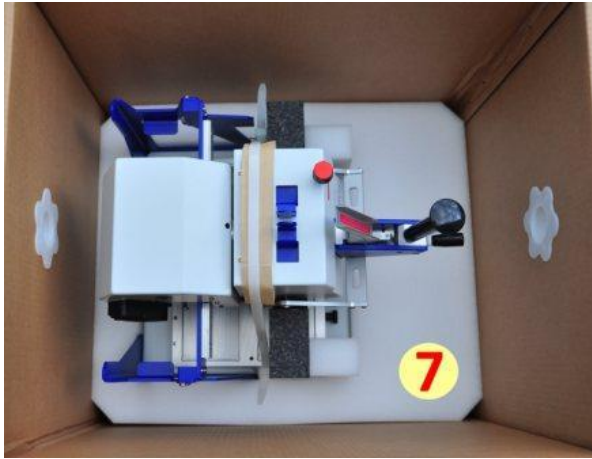

4. Insert Liquidator from above into the shipping box and place it in the lower foam piece (7). The base plate of Liquidator fits exactly into the opening of foam piece (7). Push the instrument down to assure that it sits evenly on the bottom.

5. Cover Liquidator with plastic bag (9) and place foam piece (8) over the device as shown.



6. Fill in the **Safety Declaration** and put it on top before closing the shipping box (6). Tape tight to ensure that the device is put under slight pressure and immobilized. Mark container as fragile.

Your Liquidator will only be serviced with complete safety declaration.

Liquidator is now ready for shipping. Please contact us for shipping modalities.

Steinbrenner Laborsysteme GmbH
In der Au 17
D-69257 Wiesenbach

mail@steinbrenner-laborsysteme.de

Tel: +49 (0) 6223 - 861247

Fax: +49 (0) 6223 - 861248